




Capsule Machine, ProFiller 1100, Change Part and Orienter Base Size 5

Description: Change parts include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Orienter Base / Capsule Locker

Related Products:


- Capsule Machine, ProFiller 1100, 50 Pin Tamper #5 (7472)


Installing Change Parts:





No Tools Required


Profiller 1100 system's patented design requires no tools or adjustments to change capsule size.


- 

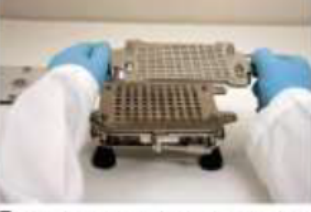
1. Unscrew 2 thumbscrews holding Orienter Tray to Base.
- 

2. Position new Orienter Tray and secure with 2 thumbscrews.
- 

3. Remove Caps Tray.
- 

4. Set Cam Lever to 3 o'clock. Remove the 3 Body Sheets.
- 

5. Lift out clear Spacer Plate and replace with new size Spacer Plate.
- 

6. Place new size Body Sheet "a" then "b", then "c" on locating pins.
- 

7. Place new size Caps Tray onto Filler. You are now ready to fill a new capsule size.

USA | Canada | Australia | International
medisca.com