



PRODUCT NO.: 7503



CHANGE PARTS, PROFILLER 1100 (Size 4)

DESCRIPTION: Change parts* for one capsule size include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Capsule Locker

ADDITIONAL INFORMATION:

*Tamper may be required. Please provide your MEDISCA Sales Representative with the FB serial # (located underneath the Filler Base) and the capsule sizes currently being filled on your current capsule machine to ensure that all pieces are included for proper capsule filling.

Related Products:

- 7457: Tamper, 50 Pin, Size 4

Installing Change Parts:

No Tools Required

Profiler 1100 system's patented design requires no tools or adjustments to change capsule size.

1. Unscrew 2 thumbscrews holding Orienter Tray to Base.

2. Position new Orienter Tray and secure with 2 thumbscrews.

3. Remove Caps Tray.

4. Set Cam Lever to 3 o'clock. Remove the 3 Body Sheets.

5. Lift out clear Spacer Plate and replace with new size Spacer Plate.

6. Place new size Body Sheet "a" then "b", then "c" on locating pins.

7. Place new size Caps Tray onto Filler. You are now ready to fill a new capsule size.

Please note that the information contained herein is for informational purposes only. The information is based on recognized compendia, manufacturers' instructions and/or other authoritative sources and texts. MEDISCA takes no responsibility for the validity or accuracy of this information. Customer MUST refer to USP, FDA, Health Canada, TGA and/or their State Board to ensure they comply with the specific regulations and ensure that appropriate procedures are in place.