

PRODUCT NO.: 7506



CHANGE PARTS, PROFILLER 1100 (Size 1)

DESCRIPTION: Change parts* for one capsule size include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Capsule Locker

ADDITIONAL INFORMATION:

*Tamper may be required. Please provide your MEDISCA Sales Representative with the FB serial # (located underneath the Filler Base) and the capsule sizes currently being filled on your current capsule machine to ensure that all pieces are included for proper capsule filling.

Related Products:

7410: Tamper, 50 Pin, Size 0-1

Installing Change Parts:





4. Set Cam Lever to 3 o'clock. Remove the 3 Body Sheets.



Orienter Tray to Base.



5. Lift out clear Spacer Plate and replace with new size Spacer Plate.



2. Position new Orienter Tray and 3. Remove Caps Tray. secure with 2 thumbscrews.



6. Place new size Body Sheet "a" then "b", then "c" on locating pins.





7. Place new size Caps Tray onto Filler. You are now ready to fill a new capsule size.

Page 1 of 1