



**PRODUCT NO.: 7508**



**CHANGE PARTS, PROFILLER 1100 (Size 00)**

**DESCRIPTION:** Change parts\* for one capsule size include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Capsule Locker

**ADDITIONAL INFORMATION:**

\*Tamper may be required. Please provide your MEDISCA Sales Representative with the FB serial # (located underneath the Filler Base) and the capsule sizes currently being filled on your current capsule machine to ensure that all pieces are included for proper capsule filling.

**Related Products:**

- 6999: Tamper, 50 Pin, Size 00

**Installing Change Parts:**

**No Tools Required**

Profiler 1100 system's patented design requires no tools or adjustments to change capsule size.

**1.** Unscrew 2 thumbscrews holding Orienter Tray to Base.

**2.** Position new Orienter Tray and secure with 2 thumbscrews.

**3.** Remove Caps Tray.

**4.** Set Cam Lever to 3 o'clock. Remove the 3 Body Sheets.

**5.** Lift out clear Spacer Plate and replace with new size Spacer Plate.

**6.** Place new size Body Sheet "a" then "b", then "c" on locating pins.

**7.** Place new size Caps Tray onto Filler. You are now ready to fill a new capsule size.