



PRODUCT NO.: 7993



CHANGE PARTS, PROFILLER 3600 (Size 3)

DESCRIPTION: Change parts* for one capsule size include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Capsule Locker


ADDITIONAL INFORMATION:

*Tamper may be required. Please provide your MEDISCA Sales Representative with the FB serial # (located on back of the Filler Base) and the capsule sizes currently being filled on your current capsule machine to ensure that all pieces are included for proper capsule filling.

Related Products:


 7582: Tamper, 60 Pin, Size 2-3

Installing Change Parts:




No Tools Required


ProFiller 3600 system's patented design requires no tools or adjustments to change capsule size.




1. Lift out Orienter Tray.




2. Slide new Orienter Tray in place.



3. Unscrew 4 thumbscrews from corners of Filler Base.




4. Remove Body Sheet Set.



5. Switch Spacer Plate.



6. Position new size Body Sheet Set on Filler Base.



7. Secure with four thumb screws.