



**PRODUCT NO.: 7995**



**CHANGE PARTS, PROFILLER 3600 (Size 1)**

**DESCRIPTION:** Change parts\* for one capsule size include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Capsule Locker

**ADDITIONAL INFORMATION:**

\*Tamper may be required. Please provide your MEDISCA Sales Representative with the FB serial # (located on back of the Filler Base) and the capsule sizes currently being filled on your current capsule machine to ensure that all pieces are included for proper capsule filling.

**Related Products:**

7984: Tamper, 60 Pin, Size 0-1

**Installing Change Parts:**

**No Tools Required**

ProFiller 3600 system's patented design requires no tools or adjustments to change capsule size.

**1. Lift out Orienter Tray.**

**2. Slide new Orienter Tray in place.**

**3. Unscrew 4 thumbscrews from corners of Filler Base.**

**4. Remove Body Sheet Set.**

**5. Switch Spacer Plate.**

**6. Position new size Body Sheet Set on Filler Base.**

**7. Secure with four thumb screws.**