



**PRODUCT NO.: 7996**



## CHANGE PARTS, PROFILLER 3600 (Size 0)

**DESCRIPTION:** Change parts\* for one capsule size include: Caps Tray / Body Sheet Set (A, B, C) / Spacer Plate / Orienter Tray / Capsule Locker


### ADDITIONAL INFORMATION:

\*Tamper may be required. Please provide your MEDISCA Sales Representative with the FB serial # (located on back of the Filler Base) and the capsule sizes currently being filled on your current capsule machine to ensure that all pieces are included for proper capsule filling.

### Related Products:


 7984: Tamper, 60 Pin, Size 0-1

### Installing Change Parts:




**No Tools  
Required**


ProFiller 3600 system's patented design requires no tools or adjustments to change capsule size.




**1. Lift out Orienter Tray.**



**2. Slide new Orienter Tray in place.**




**3. Unscrew 4 thumbscrews from corners of Filler Base.**




**4. Remove Body Sheet Set.**



**5. Switch Spacer Plate.**



**6. Position new size Body Sheet Set on Filler Base.**



**7. Secure with four thumb screws.**